

Polarization Guidelines

More than 30 minutes disconnected from a polarizing device requires **6 hours** of polarization.

Less than 30 minutes but greater than 15 minutes disconnected from a polarizing device requires **6 minutes** of polarization.

Less than 15 minutes but greater than 5 minutes disconnected from a polarizing device requires **4 minutes** of polarization.

Less than 5 minutes disconnected from a polarizing device requires **2 minutes** of polarization.

Measurement and Control Products for Science and Industry

19 Thomas, Irvine, California 92618 USA

Phone: 949.829.5555 Toll-Free: 800.288.2833

Fax: 949.829.5560 E-Mail: sales@broadleyjames.com

