

How Should A pH Electrode Be Installed?

All pH electrodes (and DO sensors) can be mounted vertically, into a vessel headplate. Electrodes and sensors can also be installed on large vessels, through a side port. For side-entry installations, the probe must be at least 15° above horizontal to insure proper operation. Remember, there is liquid inside all pH electrodes and DO sensors. This slight 15° incline is enough to prevent a bubble from lodging itself at the tip of the sensor, preventing the sensor from operating properly.

The contents of this publication are presented for information purposes only, and while every effort has been made to ensure their accuracy, they are not to be construed as warranties or guarantees, express or implied, regarding products or services described herein or in their use or applicability. All sales are governed by our terms and conditions, which are available on request. We reserve the right to modify or improve the design or specification of such products at any time.

© 2022 Broadley-James Corporation. All rights reserved. Visit www.broadley-james.com/trademarks for trademark information.

TMP-BF-102101



North America and Pacific

Email: sales@broadleyjames.com
Web: www.broadleyjames.com

19 Thomas, Irvine CA, 92618 USA

Phone: 949-829-5555
Tollfree: 800-288-2833
Fax: 949-829-5560

United Kingdom and EU Countries

Email: sales@broadleyjames.co.uk Phone: +44 (0)1525 862518
Web: www.broadleyjames.eu Fax: +44 (0)1525 862811

Wrest Park, Silsoe Beds MK45 4HS, UK